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Für MyGo Pro User **Advanced Settings:** Integration time (s) 0.5
Acquisitions per Cycle: 10

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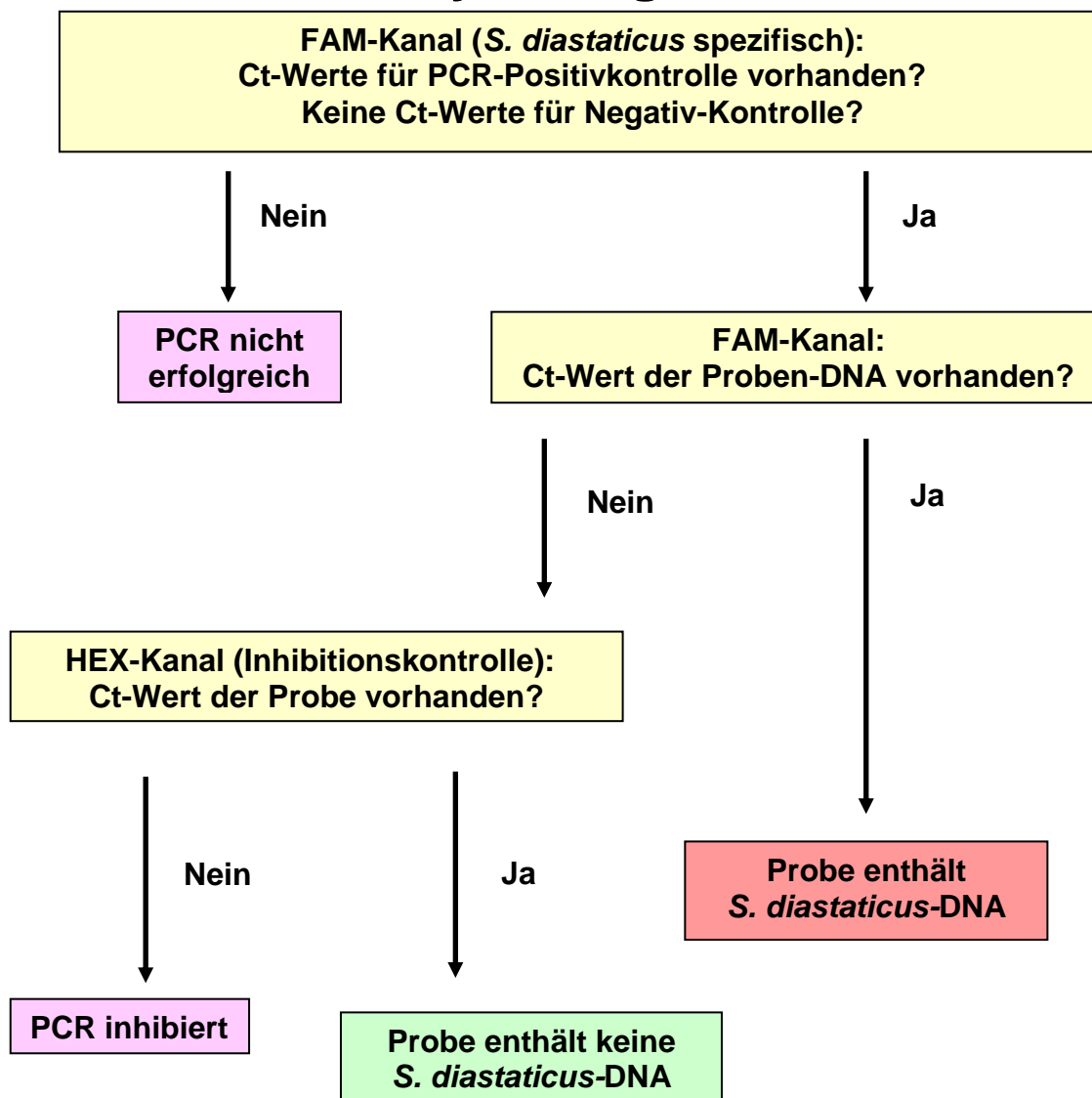
(Version 02/18)

1. Intended use

Detection of *Saccharomyces diastaticus* in wine and grape must by real-time PCR.

2. Test principle

The real-time PCR is based on hot-start-PCR and sequence-specific dual labelled probes (TaqMan[®]), which, when accurately hybridised, emit a measurable fluorescent signal of a defined wavelength in the extension phase. The increase of signal is continuously measured in a real-time PCR detection instrument.

The kit contains a specific system for the detection of *S. diastaticus*. The system emits a maximum fluorescent signal in the FAM/DQ channel. To avoid false negative PCR-results an Inhibition-Control is amplified together in the same reaction vessel with the specific sequence (HEX/DQ channel). The system contains dUTP.

Optional: Use of Uracil-N-Glycosylase will eliminate any contamination with Uracil containing amplicons from former PCRs (the enzyme is not part of this kit). **The tubes contain Lyticase.**

3. Kit contents

The kit contains sufficient reagents for 48 reactions.

1 x Premix QTPYSD	white cap
48 x QTPYSD mix (freeze-dried, incl. Lyticase and IC-DNA)	tube strips
1 x ddH ₂ O	colourless cap
1 x Control-DNA (freeze-dried)	yellow cap

4. Storage conditions

The Control-DNA is freeze-dried and has to be solved in ddH₂O prior to use (see 6.1).

Do **not** freeze the lyophilized QTPYSD mix and lyophilized Control-DNA.

All PCR reagents, except the Premix should be stored at 4 – 8 °C (39 – 46 °F).

Keep Premix for storage at - 20 °C (- 4 °F). Avoid loss of sensitivity by repeating freezing and thawing more than 3 times. For irregular use aliquot the Premix.

The QTPYSD mix contains the fluorescent labelled probes and should be handled light protected.

All reagents are stable for 6 months, if they are stored correctly

5. Materials required, but not provided

5.1. Instruments

Real-time PCR machine for low profile tubes clear

Centrifuge for 1.5 – 2.0 mL vials

Centrifuge for strips

Pipettes

“Vortex”

5.2. Reagents and plastic ware

sterile ddH₂O

sterile filter tips

optional: Uracil N-Glycosylase (0.01 U/μL added to the PCR reaction mix)

6. PCR

6.1. PCR-setup

When using the kit for the first time, the freeze-dried Control-DNA has to be shortly centrifuged and carefully resolved:

- add 55 μL sterile ddH₂O to the freeze-dried Control-DNA
- after 15 minutes mix well

Before every use thoroughly mix all PCR components and centrifuge briefly.

PCR-reaction per sample:

PCR-components	amount (μL)
Premix QTPYSD	15,0
Sample-DNA	5,0*
Total volume	20,0

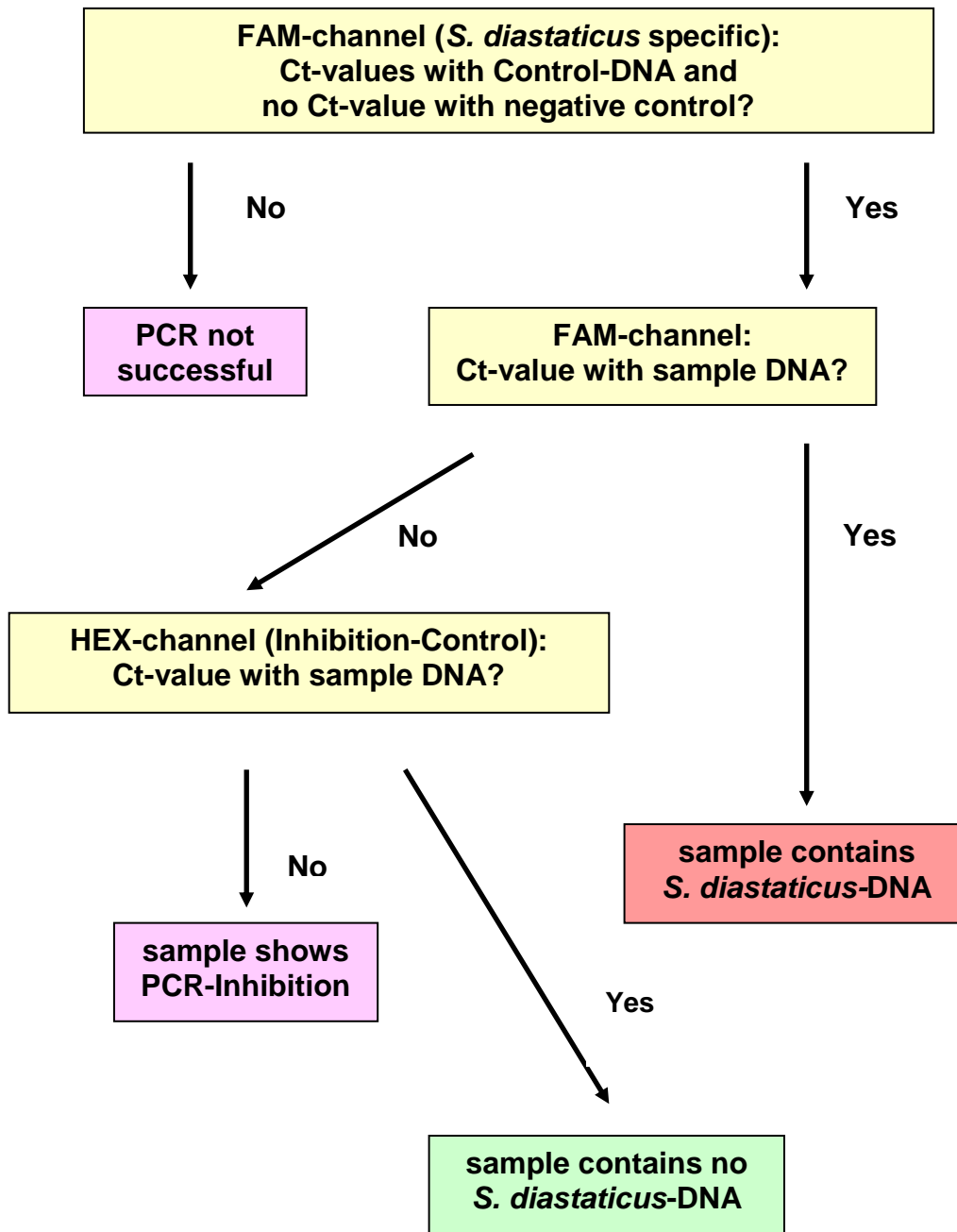
* if using the Simplex Easy DNA Kit: add 2.5 μL DNA and 2.5 μL sterile ddH₂O

1. Pipette 15 μL of the Premix into each PCR-tube, making sure that, prior to the first filling, the tip has been moistened.
2. Add 5 μL sample DNA, add 5 μL of the Control-DNA for the PCR positive control, add 5 μL of the extraction control and 5 μL of ddH₂O for the negative control* reaction. Use a fresh tip with each DNA filling.
3. Close the tubes immediately and centrifuge them shortly.
4. Place the tubes in the PCR-machine and start run.

Very important: * Please fill up the negative control with 5 μL ddH₂O to avoid unspecific amplification.

Work swiftly to avoid warming up and keep away from light.

analysis flowchart

The polymerase-chain reaction (PCR) is protected by patents and requires a licence from Hoffmann-LaRoche Inc.. The provided product does not authorise the purchaser for the commercial use of this method. GEN-IAL makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, GEN-IAL will provide a replacement product. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. GEN-IAL shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.